

Ultrospec 2000 UV/Visible Spectrophotometer User Manual

English

Deutsch

Français

Español


Italiano





Pharmacia Biotech

Keyboard

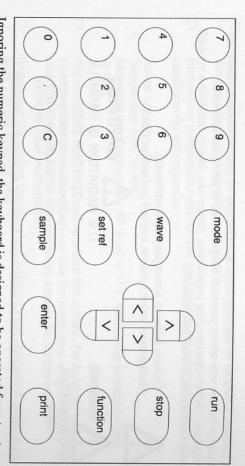


ble when the function or mode key is pressed; these options are described later. make. In addition, a range of instrument utilities and instrument modes is accessibottom and from left to right. Basic absorbance measurements are very easy to Ignoring the numeric keypad, the keyboard is designed to be operated from top to

- mode Select measurement mode required using < > \ \ \ \ keys. Press enter
- wave Select wavelength either from the numeric keypad or by using S keys. Press (enter)
- set ref units and 100 % T, respectively, on a reference solution. Set absorbance and transmittance readouts to 0.000 absorbance
- sample Select sample position in cell changer either from the numeric keypad or by using < > keys. Press (enter
- enter Select options on the display. (C) Clears a numeric entry.
- run Start measurements when operating from within the measurement modes
- stop Ends current programme
- function) Access to Set-up and Parameters.
- print Output information on the display panel to a parallel printer
- English

Display Panel

cell position number 1 in the light path. values are shown, and represent a wavelength of 360nm in absorbance mode, with The display panel has 20 characters divided up as indicated. Default start-up

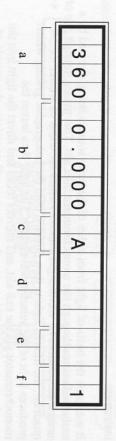

Wavelength

a)

- Display of current wavelength in nm
- 9 **Numeric Result**
- Unit of Measurement Display of numeric result unit (A is Absorbance Units; %T is Display of absorbance, % T, concentration, absorbance ratio result.

0

0 **Mode of Operation**

% Transmission; C is concentration units).

Display of instrument enhancement mode:

RAT is absorbance ratio mode; SCN is scan to chart recorder KNR is display of absorbance with time; PRO is protein concentration mode; DNA, RNA, OLI are DNA, RNA, oligonucleotide quantification modes; FC and SC are factor and standard concentration modes; KND is display of change of absorbance with time:

- e **Temperature Indicator**
- Display of temperature status if a peltier heating accessory is in use.
- °C is at temperature; LO is under temperature or heating up. HI is over temperature or cooling down;
- Sample Number

5 Display of sample number.

OPERATION

Introduction

Your UV/Visible spectrophotometer is a simple to use, microprocessor controlled

then amplified and displayed. interest and a defocussing lens to a solid state detector unit. The resulting signal is sample compartment. This passes through your cell containing the sample of monochromator via the exit slit, and mirrors focus and direct the light into the produces light of the selected wavelength. The light then leaves the quadrant; the filtered light is then directed onto the holographic grating which several (dependent on wavelength selected) filters mounted on a filter motorised mirror through the monochromator inlet slit. This passes through one of It works on the basis of light from either of the lamp sources being directed by a

Your spectrophotometer:

- measures standard absorbance, concentration and transmittance
- nucleic acid solutions. and purity checking, as well as for protein contamination measurement in has stored parameters for DNA, RNA and oligonucleotide quantification
- measures absorbance ratios
- data or output the absorbance time plot to a chart recorder (synchronised). measures change of absorbance with time and can either print the numerical
- outputs a wavelength scan to a chart recorder (asynchronous).
- of results. can be connected to a standard centronics parallel printer for output
- can be linked to a PC via its serial interface and used with the environment. SWIFT range of software application programs operating in the Windows

A range of accessories further enhances the capability of the instrument.

Measurement Modes

mode

right in the matrix shown below. Each mode is described subsequently. continuous loop. After pressing mode, use the arrow keys to move up/down/left/ The mode key gives access to the 6 operational modes; these are linked in a

Select from the table below with

mode

Concentration \$

FACTOR CONCENTRATION

STANDARD CONC

enter

Absorbance

1

Nucleic Acids

PROTEIN IMPURITY DNA OLIGO

ABS RATIO

RATE - RAW DATA RATE - DELTA DATA

SCAN TO RECORDER

Reaction Rate

1

Scan

\$

Transmission

















mode

enter to choose required measurement mode.

can return to the previous step of operation By pressing the mode button within a measurement mode the user

A flashing cursor indicates an option which is currently selected. The instrument can be switched off at any stage of operation.

Absorbance

absorbance measurements on samples, and it measures the amount of light that has been passed through a sample relative to a blank (this can be air).

Absorbance mode is the default after power on. It is used to perform simple

mode

set ref

Z S

wavelengths in the mode selected. During a standard operating to set absorbance to 0.000 AU on a reference solution at all into cell holder, by the message set reference. This is displayed procedure, the user is prompted to insert a cell containing reference

to start making measurements; sample number (and cell position) are automatically incremented after measurement.

until the reference is inserted

and set ref is pressed.

stop

to stop making measurements or return to Absorbance mode.

sample V

increments sample number (and cell position if cell changer is fitted).

if a single cell holder or sipper is installed. which is supplied with the instrument. There are different self explanatory displays The above and following sequences of instructions apply to the 6 cell changer

and can be put in sequentially if a 6 cell changer is fitted

enter

enter

(sample) with 2 samples

cell position of the <u>next</u> sample, since it is now in the light path. The advantage of using run is that if a printer is connected and RUN PRINT in Setup selected, then

there is automatic print out of the absorbance details of the sample just measured.

sample absorbance is measured and the cell changer rotated sequentially to the next

absorbance and cell position of the sample in the light path. If run is pressed the and cell position 2 selected, the display will show the wavelength used and the reference on it, remove the blank, and insert the sample(s). If sample is pressed The procedure is to select the appropriate wavelength, insert blank and set

position; the display therefore shows the wavelength used, and the absorbance and

| sample | | set ref | wave | mode |
|----------------|------------------------|--------------|---------------|-------------|
| cell | 546 | 546 | Wave | 360 |
| 11 | 546 1.200 A | 546 0.000 A | Wavelength = | 360 0.000 A |
| 2 | _ | | 546 | _ |
| cell position? | Absorbance of sample 1 | remove blank | Set λ | |

Using - R with 2 samples

546 1.412 A

2

Absorbance of sample 2

enter


360 0.000 A

| Absorbance of | 2 | 546 1.412 A | |
|---------------|-----|--------------|------|
| Absorbance of | | 546 1.200 A | |
| remove blank | _ | 546 0.000 A | tref |
| Set λ | 546 | Wavelength = | lve |
| | | | |

set

wa

enter enter

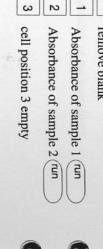


546 0.000A



00

Scan

chart recorder. Ensure chart recorder is connected, as scan data is not visible without this. The Pharmacia Biotech REC 101 is recommended An absorption spectrum can be obtained with your instrument by output to

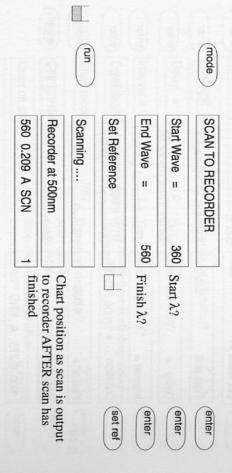


asynchronous (the chart paper drive is controlled on/off Output to the Pharmacia Biotech REC 101 chart recorder is on zero, set to internal. deflection (output is 100 mV for 1.000 Abs unit), recorder on, zero suppress Suggested chart recorder settings are 10mm/second, 200 mV full scale automatically). Chart recorder cable (80-2105-95) is required.



software module is available for use on PC. If you require post run data manipulation routines, the SWIFT-SCAN

Transmission

with that which has passed through the reference. This is displayed as a specified wavelength that has passed through the sample and compares it In transmittance mode the instrument measures the amount of light at the very high absorbances (low transmittances) transmittance mode is rarely used experimentally except for samples having its transmittance at any given wavelength is not linear, and hence percentage. The relationship between the concentration of the sample and

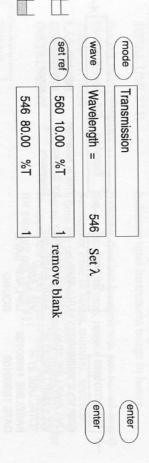


English

Messages

self explanatory, are either as follows or relate to a fault condition which requires a reason the calibration is not completed satisfactorily, the messages displayed, if not The spectrophotometer goes through a multi step calibration sequence. If for any al supplier. Other messages are displayed when an


| Display | Error Messages | service engineer from your loca accessory is installed. |
|---------|----------------|--|
| Po | | your loca |

ssible Causes

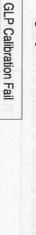

Remedy

Refer to Appendix GLP Diagnostic Tests

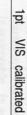

Visible region only

Check and replace UV Lamp if necessary

UV region only calibrated

1pt

V

calibrated

Check and replace visible lamp if necessary

n

Abs Non - Linear (GLP on)

Too much light

Dirty Filters compartment Too much light in sample Misalignment of filter quadrant temperature

Contact service engineer

working temperature Let instrument reach

Instrument not at working

is in place (if relevant) ensure baseplate plug Close lid properly and

Not enough light getting Check light beam is not blocked and cell compartment empty

If using a sipper: remove

to detector

Beam Blocked

PRINT OUT is ON, printer is RUN PRINT is ON, GLP off line or out of paper

Printer Error

No active printer

No Printer

Check printer state

Connect or switch on printer

or deselect RUN PRINT and GLP PRINT OUT



Temperature failed

Call Service

T

Heated accessory

unplugged

Plug in and/or press

sample

sample

error code

and advise diagnostic

Contact service engineer

Instrument warming up

Waiting for GLP

GLP Diagnostic Tests Refer to Appendix

< Min A

(<-3.0AU)

Abs value out of range

Close lid properly baseplate plug is fitted much light. Ensure Detector seeing too

> Max A

(>+3.0AU) Abs value out of range

trated. Something Replace the visible blocking light path Sample too concen-

Vis mirror fail!

enough energy during calibration Detector did not see

Accessory Messages

After installation of accessory, press

4 cell changer



(sample)

sample

Single Cell Holder

cell changer

Thermostat holder

Electric Cel

4 position cell changer

All 6 position cell changers

except ... All single cell holders,

Electric Cell Holder

Peltier Cell Holder

English

Issue 03 - 10/96