T-100 Turbidimeter

Getting Started

This instrument operates on the nephelometric principle of turbidity measurement and is designed to meet the criteria specified in ISO 7027 and DIN 27027 standards.

Using a Phillips screwdriver, remove the screws from the battery cover and install (4) AAA batteries (included). Tighten the screws to retain a waterproof seal when completed.

Keypad Functions

	Powers on and shuts off the meter. The meter automatically shuts off 20 minutes after last key press.
	Initiates the calibration mode. Press during calibration mode to abort the calibration and enter Standby "STBY" mode.
	When in measurement mode, the READ/ENTER key is used to perform a measurement. Single Measurements: When the key is pressed and released immediately (less than 0.3 seconds), the display will blink "--Rd--" for 10 times and then display the measured value. Continuous Measurements: If the READ/ENTER key is pressed and held, the instrument will perform a continuous measurement during which the display is updated every 2 seconds. This can be used for indexing vials. (This function is not available in calibration mode). After the READ/ENTER key is released, the instrument will automatically perform a single measurement.
	During calibration mode, these keys are used to select the calibration point in an incremental/decremental manner. Note: This meter uses an auto incrementing feature, i.e. after the successful calibration of one point it automatically selects the next calibration point, or automatically exits the calibration mode after the fourth calibration point. See also "Restoring Factory Calibration".

Turbidity Calibration

The T-100 was calibrated and tested prior to leaving the factory. The T-100 kit includes (4) primary standards (0.02, 20.0, 100 and 800 NTU) to verify the meter calibration or to recalibrate the meter as needed. Before measuring samples for the first time, perform a meter calibration with the primary standards. Four standard vials are included and labeled for each standard level. 800 NTU (CAL 1), 100 NTU (CAL 2), 20.0 NTU (CAL 3), 0.02 NTU (CAL 4).

Place the light shield cover over the vial to keep stray light from changing the calibration and sample measurement readings. In many environments, the light shield is not required as the vial cap is able to keep stray light from entering the sample well.

Indexing is not required due to the high quality of the glass vials. Simply align the index mark (arrow) on the vial with the index mark (arrow) on the meter.

Note: Do not shake the calibration standards as air entrapment will introduce error. Do not allow the calibration standards to freeze as this will cause permanent damage to the standards, making them inaccurate.

Calibration Procedure

Preparing the Calibration Standard Vials


Ensure there are no scratches or marks on the glass vial. If there are marks that could affect the turbidity measurement, select another vial. Before using the standards for the first time, rinse the inside of the vial with a small amount of the same NTU standard that will be used to fill that vial.

1. Gently pour about 5 mL from the 60 mL stock standard into the vial.
2. Cap and swirl the vial so the inside of the vial is fully rinsed with the standard.
3. Pour the standard out of the vial and discard it.
4. Repeat steps 1 through 3.
5. Gently fill the vial using the 60 mL stock standard above the fill line on the vial and cap the vial.
6. Repeat this rinse and filling procedure for the other three vials.

Note: The vial rinsing is critical for precision with the lower NTU standards.

Use these calibration standard vials daily as a check, and also periodically for calibration as needed. When the standard reading of 20.0, 100, or 800 NTU drifts more than 10 % of the standard's NTU value, the solution should be replaced with another aliquot from the plastic bottle following the steps listed above. **Note:** 0.02 NTU standard read back value should be <0.1 NTU.

Calibration Sequence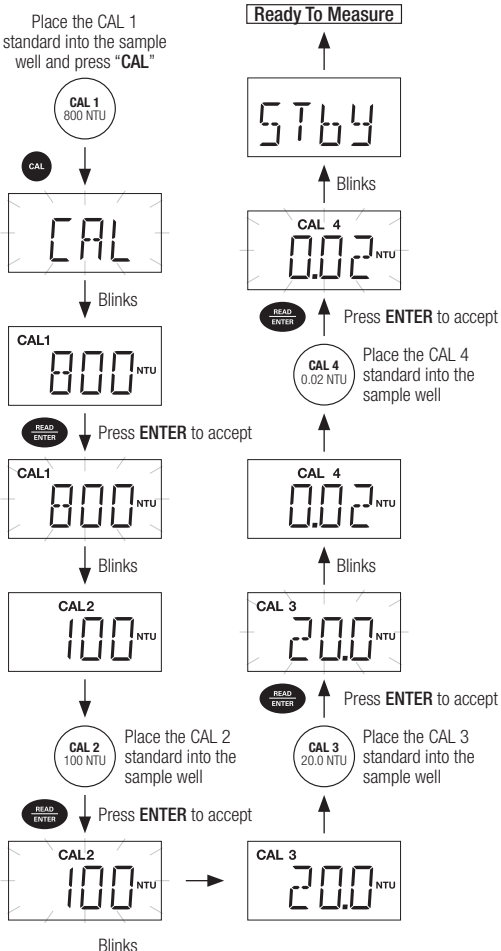


Restoring Factory Calibration

1. With the meter off, press and hold key.
2. Power on the meter using **ON/OFF**. Release key when "URST" (User-Reset) appears in the display.
3. The display shows "URST" for about 2 seconds and then shows "No".
4. Use or key to choose "YES" or "No".
YES = To restore meter back to factory calibration values
No = To retain last calibrated values
5. Press **READ/ENTER** key to confirm. When "YES" is selected, the meter performs a reset, and will blink "--Rd--" to indicate that it is reading in measurement mode.

Turbidity Measurement

Insert vial containing sample or calibration standards before powering on.

Alternatively, from "STBY" mode, press **READ/ENTER**.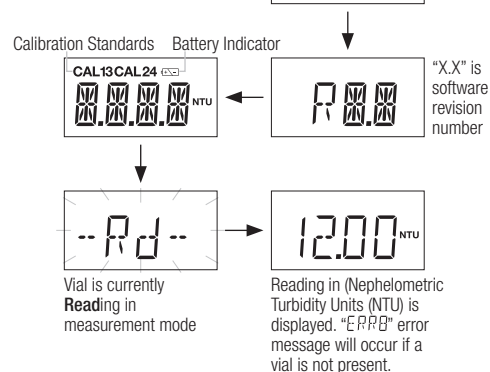


Turbidity Measurement

Good techniques are important for accurate turbidity measurements. For best results, use clean sample vials, properly position the vial in the sample well, cover the vial with the light shield cover, use proper handling of standards, use supplied cloth to keep vials clean and scratch free, and perform meter calibration as needed.


Preparation of Sample Vial

1. Obtain a clean, dry sample vial.
2. Handle the sample vial only by the cap to avoid fingerprints on the glass.
3. Rinse the vial with approximately 10 mL of the sample, capping the vial with the black screw cap and gently inverting it several times. Discard the used sample and repeat.
4. Fill the rinsed vial with the sample up to the mark indicated in the vial. Cap the vial with the supplied black screw cap.
5. Wipe the vial with the soft, lint-free cloth supplied. Ensure that the outside of the vial is dry, clean and free from smudges.
6. Apply a thin film of silicone oil (supplied) on the sample vial (see image below).
7. Wipe with a soft cloth to obtain an even distribution over the entire vial's surface. The purpose of oiling the vial is to fill small scratches and to mask the imperfection in the glass. Do not apply large quantity of oil as this may collect dirt and dust.
8. The sample vial is now ready to be inserted into the sample well for measurement.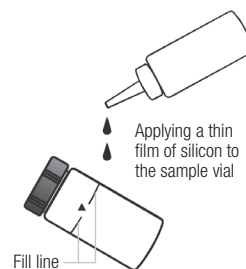

Measurement Procedure

1. Turn meter off, and place on a flat, level surface.
2. Place the sample vial inside the sample well and align the vial's index mark with the meter's index mark.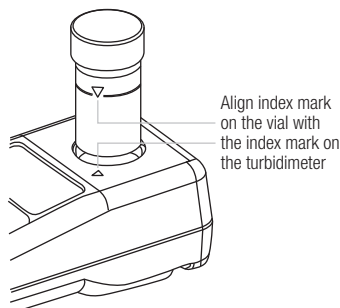


3. Push the vial down until it is seated.
4. Place the light shield cover over the vial that is inserted into the sample well.
5. Turn on the meter by pressing the **ON/OFF** key. After the power-up sequence, the meter goes to measurement mode and the display blinks "--Rd--". The measured reading appears in the display.
* When meter is in "STBY" mode after calibration it is ready for measurement. Press the **READ/ENTER** key. The display blinks "--Rd--" until the measured reading appears.

Notes:

- NEVER pour liquid directly into the sample well of the instrument. Always use a vial. The instrument will only accurately measure the turbidity of a sample when vials sealed with the black caps are used. The black cap serves as both seal and a primary light shield.
- NEVER attempt to clean the sample well. The optics may be damaged.
- For battery conservation, the instrument automatically powers off 20 minutes after the last key pressed.

Measurement Sequence

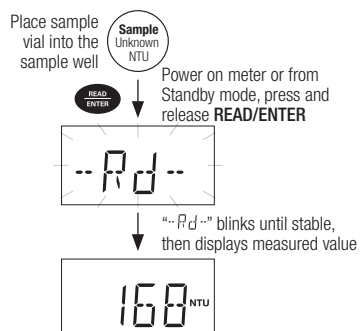Troubleshooting Guide

The T-100 turbidimeter routinely performs self-diagnostics, and will automatically generate messages to provide you with specific diagnostic information.

Message	Description	Corrective Actions
ERR1	Calibration Error. The meter is unable to recognize the 800 NTU Calibration Standard.	Check or replace the 800 NTU calibration standard.
ERR2	Calibration Error. The meter is unable to recognize the 100 NTU Calibration Standard.	Check or replace the 100 NTU calibration standard.
ERR3	Calibration Error. The meter is unable to recognize the 20.0 NTU Calibration Standard.	Check or replace the 20.0 NTU calibration standard.
ERR4	Calibration Error. The meter is unable to recognize the 0.02 NTU Calibration Standard.	Check or replace the 0.02 NTU calibration standard.
ERR5 ERR6 ERR7	General Calibration Error. There is not sufficient signal to achieve the appropriate resolution.	Perform recalibration with all (4) standards.
ERR8	Excessive stray light detected.	Ensure the vial is properly seated in the sample well. Use light shield cover.
ERR9	Lamp Failure.	Return unit.
OR	"Over Range"; value is above maximum measurement range.	Dilute sample.
	Low battery indication. The batteries need to be replaced.	Replace batteries.

Routine Maintenance and Care

The supplied carrying case is optimal for protecting the instrument. If you do not plan on leaving the instrument in the supplied carrying case, when not in use, ensure that the instrument has been turned off and that a clean sample vial fitted with a black cap has been placed in the sample well. This will ensure that a minimal amount of dust and/or debris will be able to settle on the optics of the instrument.

Proper measurement of the turbidity of a sample requires the use of a vial that is free of marks, smudges, scratches and bacterial growth. Sample vials must be handled with absolute care to avoid contamination or damage, which might change the optical characteristics of the glass. Scratches, fingerprints, and water droplets on the sample vial or inside the sample well can cause stray light interference leading to inaccurate readings.

Cleaning the vial is accomplished by washing the interior and exterior of the vial in a detergent solution. Once cleaned, the vial should be rinsed thoroughly 8 to 10 times with clean distilled water to eliminate the possibility of detergent buildup and streaking.

Vials can also be acid washed periodically and coated with a special silicone oil to fill small scratches and mask the imperfections in the glass. Since the silicone oil required for this application should have the same refractive characteristics as glass, it is recommended that the oil be obtained from us. Care should be taken not to apply excessive oil that could attract dirt or contaminate the sample well of the meter. Once the oil has been applied to the vial, the excess oil should be removed with a lint-free cloth. The result should be a sample vial surface with a dry appearance, but with all imperfections filled with oil.

Sample vials should always be handled from the top or by the cap to avoid fingerprints or smudges. After a vial has been filled with a sample and capped, the outside surface should be wiped with a clean, lint-free absorbent cloth until it is dry. Cleaned and dried vials should be stored with the black caps on. The vials can be stored in the carrying case. During normal operation you may use any typical glass cleaner along with a lint free cloth or tissue (Kimwipes®), to clean the outside of the vials.

Condensation may appear on the vial when your sample is very cold and the relative air humidity is high. When this happens, the turbidity that you read may be higher than the actual turbidity due to the light scattered by the condensate on the vial. To alleviate this problem, coat the vial with an anti-fogging agent, or by running warm water over the vial for a short period of time to warm the sample prior to measurement.

Intended Use, Maintenance & Precautions

This instrument is designed for turbidity measurement of water-based solutions. For routine maintenance dust or wipe the display using a damp cloth. If necessary, warm water or a mild water based detergent can be used. Immediately remove any spilled substance from contact with the meter using the proper cleaning procedure for the type of spill.

- Do not use this equipment in potentially explosive atmospheres.
- Ensure that no liquid enters the instrument.
- Do not use any aggressive cleaning chemicals (solvents or similar agents).
- There are no user serviceable parts inside. Attempts to service internal parts may void the warranty.
- Not intended for medical applications or patient use.
- **WARNING:** No modification of this equipment is allowed.

Instrument Operating Conditions

Operating Ambient Temp.	0 to 50 °C
Operating Relative Humidity	0 to 90 %, non-condensing at 30 °C
Storage Temp.	-20 to +60 °C (instrument only)
Storage Relative Humidity	0 to 90 %, non-condensing at 30 °C
Pollution	Degree 2
Overvoltage	Category II
Weight	200 g
Size (L x W x H)	15.5 x 6.8 x 4.6 cm
Battery Requirement	4 x AAA (LR03) 1.5 V batteries
Vibration	Shipping/handling per ISTA #1A
Shock	Drop test in packaging per ISTA #1A
Certifications/Compliance	FCC part 15 class A, TUV 3 in 1, CE, RoHS

Ordering Information

35635-00	Waterproof T-100 handheld turbidimeter with set of 4 calibration standards (800, 100, 20.0 & 0.02 NTU), sample vials, lint-free cloth, silicone oil, batteries, and hard carrying case
35635-05	Waterproof T-100 handheld turbidimeter, sample vials, lint-free cloth, silicone oil, batteries, and hard carrying case
35635-52	Calibration kit set – standard solutions (0.02, 20.0, 100, 800 NTU) 60 mL each and (4) vials
35653-55	Pack of 3 sample vials
00159II	Silicone oil (10 mL)
00131SH / 32X366401	Replacement light shield cover
67X357201	Replacement cloth
09376-00	AAA batteries 12/pack

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