ALCATEL 2008A Users Instruction Manual

- POMPE PRIMAIRE MECANIQUE
- MECHANICAL VACUUM PUMP
- MECHANISCHE FEINVAKUUMPUMPEN

I. CARACTERISTIQUES TECHNIQUES**I. TECHNISCHE DATEN**

3

I. TECHNICAL CHARACTERISTICS

3

11. UTILISATION

4

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II. OPERATION

2.1. Description

Figure 01 shows a diagrammatic cross-section and a side-view of the pump unit.

- A. Pumping section : mechanical, 2-vane, 2-stage pump
- B. Transmission : coupling
- C. Drive motor
- D. Motor switch
- E. Inlet
- F. Oil filter plug
- G. Oil level sight glass
- H. Oil drain plug
- I. Gas ballast valve
- K. Outlet

2.2. First starting

Remove the protective caps from the inlet and outlet pipes.

Connect the motor, after checking the supply voltage.

Three-phase models

Connect the terminals in the terminal box as shown in the diagram reproduced inside the lid, for 220 V or 380 V (50 Hz), 220 V or 440 V (60 Hz), as required.

Momentarily apply power and check the direction of rotation of the motor. The end of the shaft must turn in the direction of the arrow on the frame.

Single-phase models

The pump is delivered with a 2 meter long cord (2 wires plus ground). Wiring is for 220 V operation (50 Hz) or 110 / 220 operation (60 Hz).

The rotation direction is set at the factory.

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Motor protection for the rated current should preferably be provided.
Unscrew the oil filler plug and fill to the middle of the sight glass. This operation should be carried out with the pump off.

It is quite normal for the level to vary when the pump is rotating.

Oil level is always checked with pump at rest!

Any of the following oils may be used :

ALCATEL VP1 oil
SHELL TALPA 30 oil
ANTAR PV 600 oil
CODITEC PV 195 A oil

- I install the pump in the system and switch it on.
- The pump can be stopped without introducing air, as it is airtight and oil migration is impossible

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2.3 Normal starting

Since the pump is airtight, no precautions need be taken on «cold starts», provided that the temperature is not below 10°C (40°F)

2.4. Inlet connection

The pump inlet port is equipped with a NW 25 flanged pipe (P/N 52594). Connection of different nipples or fittings is secured by quick connect clamp. (see Alcatel catalogues : Vacuum Fittings, High vacuum Valves).

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If the pump has absorbed a considerable amount of condensable vapours, the condensate may become emulsified with the pump oil. In this case, it is merely necessary to close the inlet and run the pump by itself with the gas ballast valve fully open for approx. 30 min., to regenerate the oil.

2.6. Exhaust of pumped gases

The exhaust port of the pump is equipped with a **NW 25** flanged pipe (P/N 52595).

To pump out the oil mist delivered when the pump is working at high pressure a hose nipple can be connected to this flanged pipe (see 2.4.)

An oil mist eliminator (P/N 52599) can also be fitted on the outlet.

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111. MAINTENANCE

3.1. Draining the oil

The quality of the pump oil is extremely important for proper operation. The vacuum pump oil must be inspected periodically for contaminants and oil level which must be maintained in the center of the sight glass. Oil level is checked with the pump not rotating because the level will vary during operation. Alcatel pumps may be turned off for visual oil level inspection, while backing diffusion pumps etc.. as they seal under vacuum when turned off. The oil consumption due to exhausting gases will depend upon frequency of pump down and operating pressure.

Alcatel Exhaust Mist Eliminators will reduce loss considerably by condensing expelled oil vapors, and returning the oil to the reservoir.

Inspection of the oil level will also indicate oil contamination. The oil should have a light amber color. If brownish in color, the oil should be changed. Another indication of oil contamination, is an increase in base pressure of the pump. Alcatel pumps are equipped with gas ballast, which can be used to regenerate oil contaminated with condensable vapors. The oil will be cloudy with a high base pressure. In this case, close the inlet and open the gas ballast valve for approximately thirty (30) minutes. The base pressure should reduce to a normal value after closing the gas ballast valve.

3.2. Flushing

If the base pressure is still high, an oil change must be performed. The oil should be drained when hot and discarded. When the oil has drained, vent the intake and run the pump for several seconds to expel oil trapped inside the pump stages. Re-charge with fresh clean oil through the fill plug, to the middle of the sight glass.

In some extreme cases, deposits have built up inside the pump so that new oil is contaminated within a few hours. In this case, the flushing procedure should be performed, as detailed herein.

Drain oil from hot pump and re-install the drain plug. Place a cloth lightly over the exhaust port to contain oil splash. Turn the pump on, with inlet open and slowly pour fresh oil into the inlet - equal to the pump's capacity. Turn off pump and drain flushing oil. Re-charge oil level to middle of sight glass by way of the oil fill plug. Alcatel oil VPI is recommended for re-charging or flushing Alcatel vacuum pumps.

The cleanliness of the oil will greatly prolong the life and performance of the pump. A blank statement for oil changes cannot be made for all applications due to the differences in operating conditions.

A general rule is the oil change every 2 to 3 months in clean applications ie, diffusion fore pump on UHV systems.

Extreme applications, such as ion implantation, chemical stripping or etching systems could require weekly maintenance changes or expensive trapping techniques.

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3.3.1. Dismantling

The only regular maintenance required by our pumps is changing the oil or flushing.

However, prolonged absorption of dust or acid-vapor-laden gases may produce clogging which can no longer be removed by flushing out with oil.

In that event the pump must be dismantled and all parts, except gaskets and seals, washed in trichloroethylene, or perchloroethylene.

ALCATEL pumps are designed for easy dismantling, and no adjustments are required on reassembly. All parts whose relative positions must remain constant are dowelled after accurate setting at the factory, and on reassembly the various parts will always be located in their correct positions by merely tightening the screws and nuts.

The only tools required are a few wrenches, which will generally form part of the standard equipment of any maintenance department.

Before dismantling the pump, drain the oil (see section 3.1.).

The first step of dismantling (fig. 02) is to remove the pumping section from its chamber.

The second step is to dismantle the pumping section (fig. 03).

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**3.3.1. Dismantling of the pumping section
(fig. 02)**

Unscrew the four handle plate fixing nuts 71211 and the fixing screw 75530 .

Pull out the handle plate 52547 together with the motor . Be careful not to damage the flat gasket 52555 . One half of the coupling assembly remains on the pump side (fan), the other one comes out on the motor shaft.

The pumps equipped with an American made motor get an adaptor flange 52589 mounted between the handle plate and the motor.

Remove the oil casing being careful not to damage the flat gasket 52548

Unscrew the inlet outlet flange fixing screw 82734 and remove the flange 52566

Remove the inlet port 52594, the filter 52565 and the outlet port 52595.

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3.3.2. Dism antling of the pumping section (fig. 02 + 03)

Before dismantling, carefully note the positions of the bearing plates and the stators. These are precision-located on the bearing plates on assembly at the factory, and must be replaced in the same positions on re-assembly in order not to damage the dowel pin holes. Note part numbers.

Remove the cover 52.538, by unscrewing the two fixing screws 75508.

Remove the valve spring 52541 and the valves 52540

Remove the pump base 52521 by inscrewing the two set screws 75514

Put the pump upside down then turn the fan by hand and unscrew the shaft key set screw 75554

Remove the fan coupling 52552 and the key 83649.

Place the pump vertically. Remove the impeller 83290 which is fixed on the shaft by means of circlip 71161. There is no need to remove the valve lever assembly 72578 from its bracket 52567.

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a) Removal of the rear bearing plate 52571

Unscrew the three nuts 71214

Unscrew two screwdrivers into the slots at 180° to each other. Turn them about their axis to release the bearing plate from its doweling pins.

Pull out the plate along its axis, being careful not to damage the shaft seal lips 83645. Remove the o-ring 79030.

The anti-suckback system is set at the factory. If dismantled, it must be reset as described under «re-assembly». However, the rear bearing plate can be removed without disturbing the setting of the sealing system.

b) Removal of the high pressure rotor 52510

Make it slide along the double bore of the stator.
Remove the vanes 83946 and the springs 53040.

c) Removal of the high pressure stator 52572

Use two screwdrivers as indicated hereabove (a)
Remove the o-ring 79030.

d) Removal of the central bearing plate 52573

Use two screwdrivers as indicated hereabove (a)
Remove the o-ring 79030

e) Removal of the low pressure rotor 52512

Remove the rotor 52512 by sliding along the double bore of the stator, being careful not to turn it in order to prevent damaging the lips of the shaft seal 83645.
Remove the vanes 53106 and their springs 53040.

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3.3.3. Cleaning of parts and replacement of gaskets

a) Cleaning

All metal parts must be cleaned with trichloroethylene, using a brush. Rags and cloths must on no account be used.

Gaskets must not come into contact with trichloroethylene.

b) Replacement of gaskets

Check that the gaskets are in good condition and replace as required.

3.3.4. Re-assembl y

Before re-assembly, dry all parts cleaned with trichloroethylene by means of compressed air. Re-assembly must be carried out in the exact reverse order of the dismantling operations, noting however the following points :

All the metallic parts and the lips of the shaft seals 83645 will be coated with lubricating pump oil, being careful to not fill the doweling pin holes.

After the pump has been re-assembled, it is important to fill up the two holes of the front bearing plate 52575 with oil.

Use a torque wrench to tighten the three nuts 71214. Maximum torque : 1 Nm.

The anti-suction valve assembly must be readjusted only if it has been completely dismantled. Proceed as follows (see figure 04).

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a) Slowly turn the pump shaft clockwise until the axes of the impeller (A), clips (B) and the seat (C1) are in line.

b) Offset the cam while holding the impeller with a finger so as to exert a slight force in (F) direction.

With the cam offset, the anti-suckback valve should be able to rise to about 1 mm (.025").

The surface of the seat (C) and the anti-suckback valve (3) must be parallel. This adjustment is made by loosening the screw (4) and rotating seat (C).

After these operations, check that the lever returns properly to its seat (C).

3.3.5 Ordering spare part

For quick delivery of spare parts always note:

- 1) The type of pump
- 2) The serial number
These particulars are given on the specification label attached to bracket 52545.
- 3) the part number as shown in the accompanying illustrations.

To simplify maintenance, we provide ready-packed kits of most commonly required parts. These kits are always available. The use of these kits is recommended, as you are then certain of having all the parts that you are likely to need. In addition, you will receive faster service without risk of error or confusion.

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Kit of gaskets (P/N 52610)

This kit contains all gaskets included in the pump and all parts which have to be replaced whenever the pump is completely dismantled.

Two-years maintenance kit (P/N 52613)

This kit comprises, in addition to the parts in the gasket kit 52610, a set of parts required for maintenance of the pump over a period of two years, under normal working conditions.

Note : The small anti-suck back valve seat mounted on lever 52578 (supplied as an assembly) is also available separately under PN 83772.

3.6. Tools

Tools required for dismantling and reassembling the pumping unit:

Open wrenches : 13 and 10 mm
Allen keys : 5 and 6 mm
Circlips pliers i.d. : 26 mm

ILLUSTRATIONS

figure 01 - pump unit.
figure 02.1 - diagram of motor pump unit, three-phase models
figure 02.2 - dismantling of motor pump unit, single-phase models
figure 02.3 - dismantling of pumping section.

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



ILLUSTRATIONS

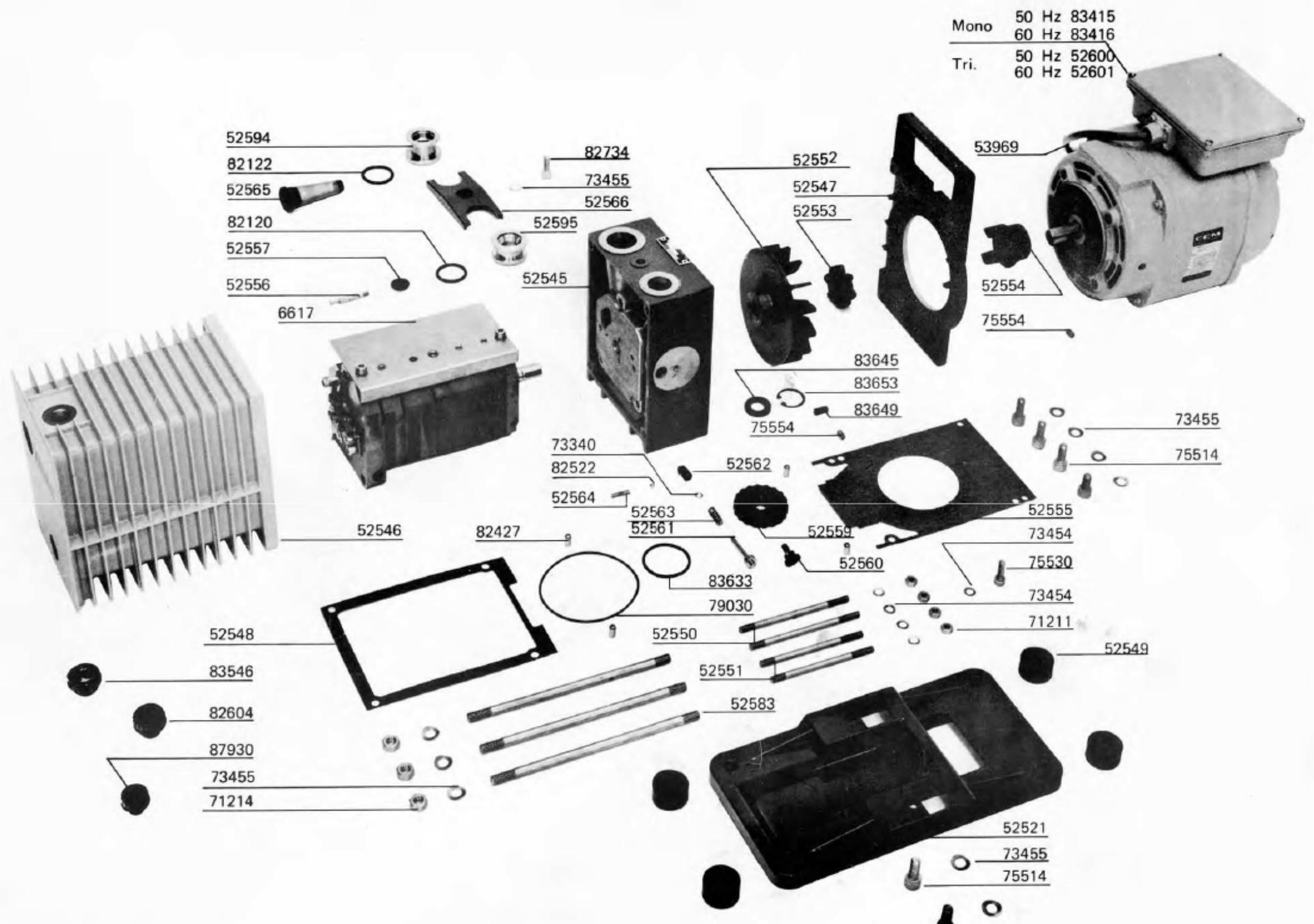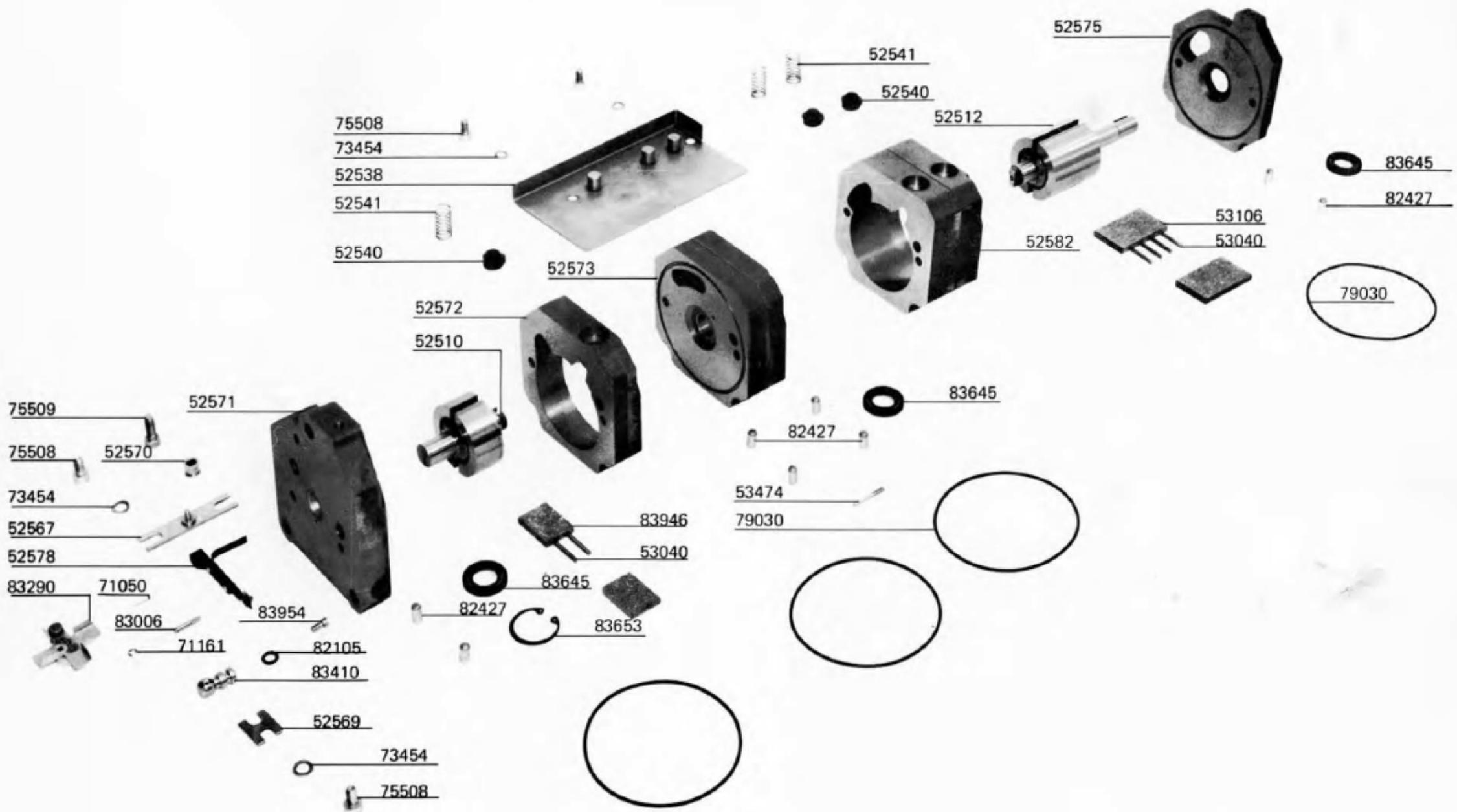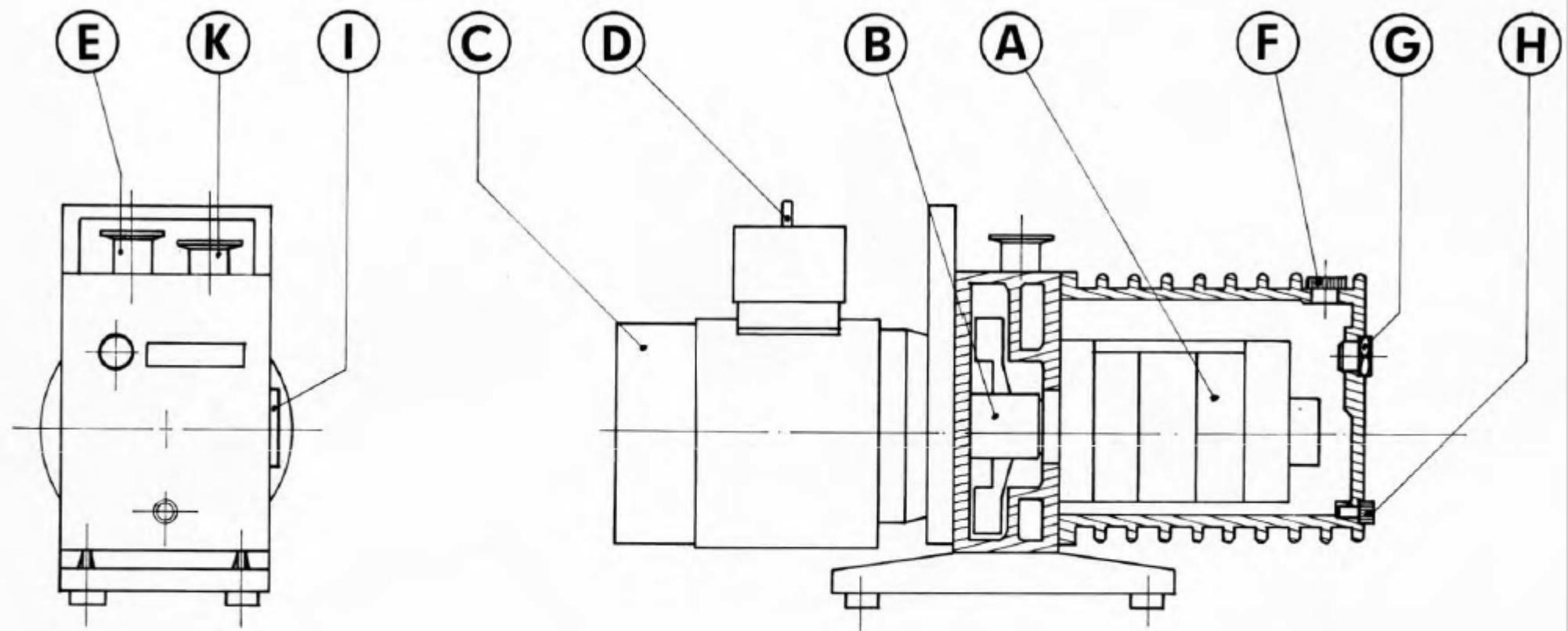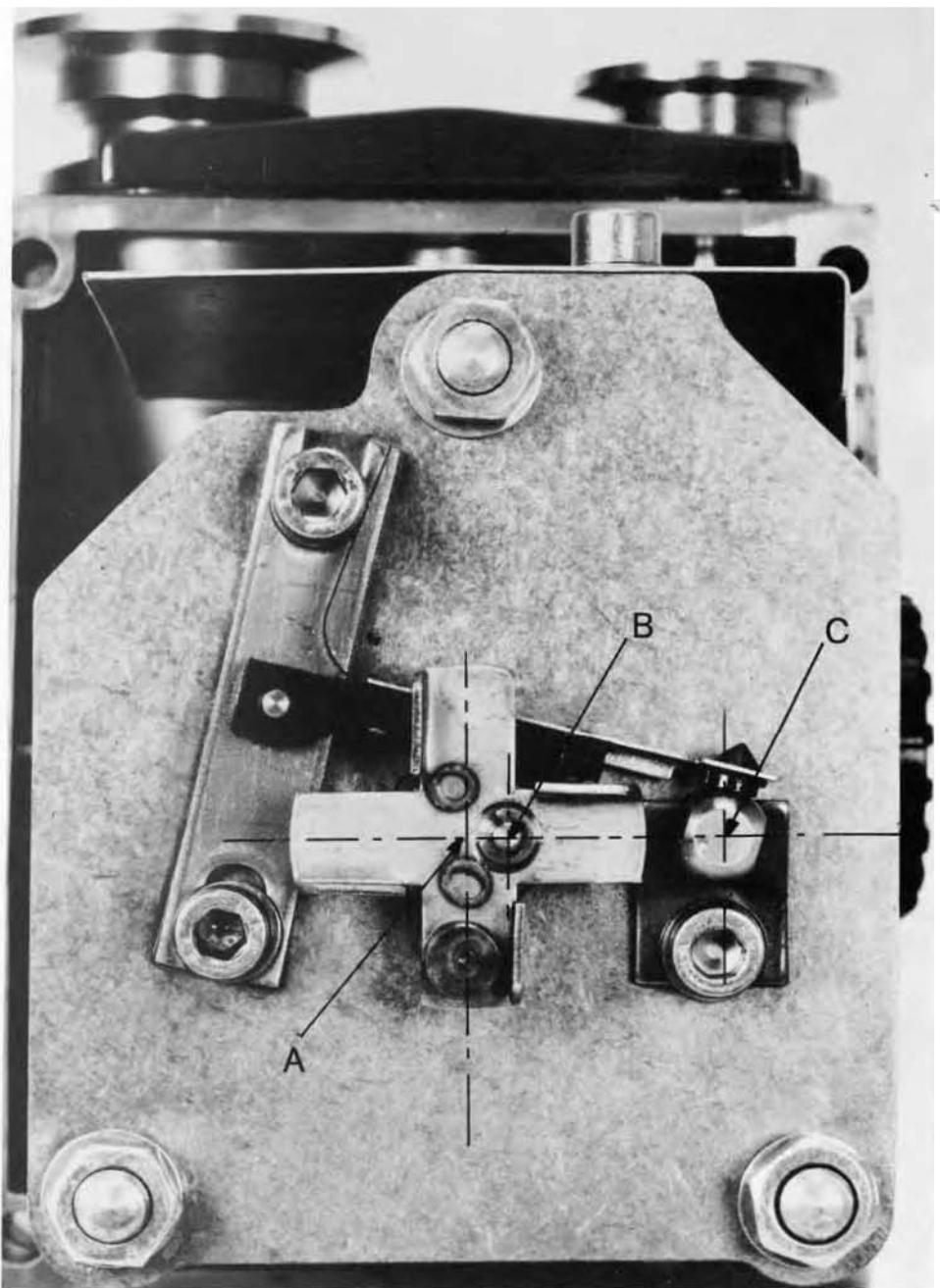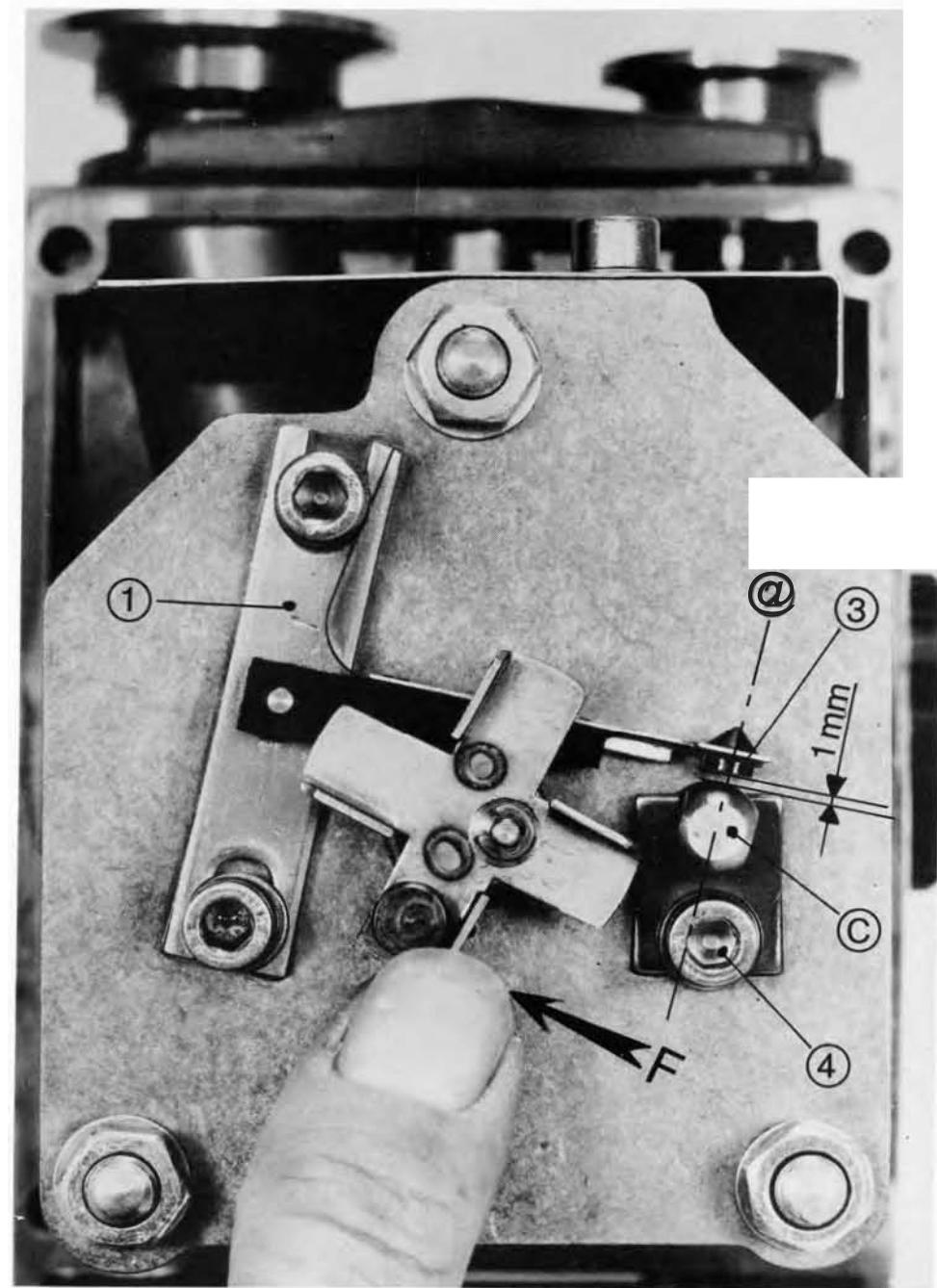
figure 01 - pump uni t.
figure 02 - dismantling of motor pump uni t.
figure 03 - dismantling of pumping section.
figure 04 - Sealing system .

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