Caframo®

Compact Digital Stirrer User Manual

Reversing Digital Stirrer User Manual

Online Warranty Registration

To register the warranty for your Caframo stirrer online, please complete the registration found at

CaframoLabSolutions.com/product-registration/

Language Index

English	1
French	17
Spanish.....	33
German.....	51

Contents

Packing List	2
Introduction	2
Safety Label Legend	2
Safety Instructions	3
General Instructions	4
Approvals	4
Installation	5
Environmental Operating Conditions	5
Auto Start.....	6
Torque Capabilities	7
BDC2002 Dimensions	8
BDC2010 Dimensions.....	9
BDC2002 Technical Data.....	10
BDC2010 Technical Data.....	11
BDC2002 General Operating Instructions.....	12
BDC2010 General Operating Instructions.....	13
Cleaning and Service	15
Warranty and Product Liability.....	16

Packing List


Compare the contents of the shipment to the items on the following list to ensure all parts are received with the stirrer. Do not discard the container and packing materials until all parts are accounted for. It is recommended that the original packing materials be kept for transportation and storage.

- Overhead Stirrer
- Power Supply with 120 volt cord set or
Power Supply with 220-240 volt cord sets (U1)
- Support Rod (including cap screw and lock washer)
- Chuck Key
- Instruction manual

U1 Cord Sets

Included with BDC2002U1, BDC2010U1.

- a. United Kingdom
- b. Europe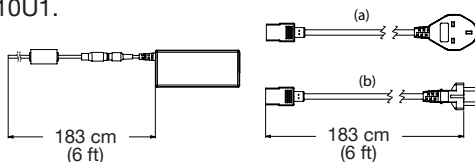

RoHS compliant & CE marked.

Introduction

This stirrer is specifically designed for professional use in the laboratory and for use in industry to stir liquids with various properties and applications.

This stirrer uses a high efficiency brushless DC motor. The speed is kept constant with continuous feedback to the motor. When torque demands are outside the stirrer's range, integrated electronic overload protection ensures no damage can occur to the electronics, motor or transmission.

The stirrer is designed for reliability and durability. There are no wear parts that require user service during the normal life of the stirrer. There are no user serviceable parts.


Safety Label Legend

CAUTION: Read operating instructions carefully and in full.

RECYCLE: Recycle in accordance with the waste electronic and electrical equipment (WEEE) directive.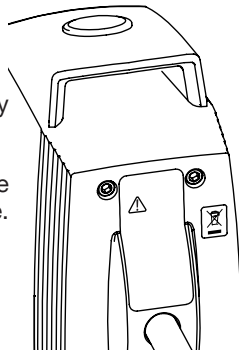

Safety Instructions

-  1. All operators must be familiar with the stirrer and read this entire manual. If the stirrer is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.
-  2. The stirrer must be securely fixed to a stable support. Mount to a Caframo stand using a Caframo heavy duty clamp. If other stands or clamps are used, the stirrer must be secured such that it will not fall if tipped.
-  3. **CAUTION:** This is not an explosion proof stirrer. Do not use in a highly flammable or explosive environment.
-  4. Spinning impellers can cause severe personal injuries. Operators must use extreme care and good judgment when mixing at any speed.
-  5. All mixing paddles and impellers must be in good condition with straight shafts. If the stirrer vibrates at high speeds check the paddle shaft for damage and repair or replace if needed.
-  6. Extreme care must be taken when mixing chemicals to ensure that no chemicals are splashed outside the mixing vessel. Care must be taken when changing to faster mixing speeds.
-  7. Ensure that the mixing impeller does not contact the containment vessel.
-  8. Do not operate while standing in water. Keep the stirrer dry and do not immerse any part, except the mixing paddle into any liquids. Protect from splashing.
-  9. Ensure that no loose clothing, jewelry, or hair can become entangled in any rotating parts. A fast spinning chuck can cause injury to operator.
-  10. Power can be interrupted to the stirrer by pressing the power button. In an emergency, disconnect the mains cord.
-  11. Ensure the power button and mains power cord of the stirrer can be accessed immediately, directly and without risk at any time, and above all, in case of fault.
-  12. The stirrer must be used with supplied power cord. The cord and grounded plug type are important safety elements. Do not use an inadequately rated cord or ungrounded plug. Always connect the mixer to a grounded outlet. Do not alter the plug.

General Instructions

Read Safety Instructions: Unpack the stirrer carefully. Ensure the stirrer is not damaged and verify the contents against the packing list. If the stirrer is damaged or the contents do not correspond with the packing list, notify your supplier immediately.

Read the manual carefully! Make sure that all users read this manual thoroughly before operating the stirrer. For further questions or inquiries, please contact:

Caframo Limited

501273 Grey Road 1
Georgian Bluffs ON N0H 2T0, Canada

Tel: 519-534-1080
Toll Free: 800-567-3556 (within USA and Canada)
Website: CaframoLabSolutions.com
E-mail: labsales@caframo.com

Approvals

Standards		
North America		CAN/CSA C22.2 No. 61010-1-12 CAN/CSA-C22.2 No. 61010-2-051-15 UL 61010-1 (3rd Edition) UL 61010-2-051 FCC Part 15, Class A ICES003
Europe, Australia, New Zealand		IEC 61010-1, IEC 61010-2-051 EMC/EMI standard IEC 61326-1

CSA	Canadian Standards Association
CE	Communité Européen (European Community)
UL	Underwriters Laboratories
FCC	Federal Communications Commission
IEC	The International Electrotechnical Commission
ICES	Interference Causing Equipment Standards
EMC	Electromagnetic Compatibility
EMI	Electromagnetic Interference

Installation

Installing the support rod

Remove the hex screw and lock washer from the support rod. Insert the support rod into the back of the stirrer so that the holes line up for the hex screw. Secure with hex screw and lock washer and tighten with hex key.

Clamp the stirrer to a Caframo stand on a stable, level surface (see Safety Instructions).

Connect stirrer to supplied Caframo power supply. Connect power supply to a grounded electrical outlet.

Impeller attachment

Insert the impeller shaft into the open chuck and adjust the impeller to the desired height. The hollow shaft allows the impeller shaft to come through the top of the stirrer if necessary. Tighten the chuck's outer ring by hand and then secure with the chuck key.

CAUTION:

 Operating a freely rotating impeller shaft through the top of the stirrer is not safe. Do not allow the impeller shaft to touch or extend through the rubber gasket while the motor is powered on. The gasket is for sealing purposes only.

Environmental Operating Conditions

The BDC2002 and BDC2010 stirrer must operate in the following conditions:

Indoors

1. Altitudes up to 2000 m (6500 ft).
2. Temperatures from 5°C to 40°C (41°F to 104°F).
3. Maximum relative humidity 80% for temperatures up to 31°C (88°F) decreasing linearly to 50% relative humidity at 40°C (104°F).
4. Pollution degree 2 in accordance with IEC 664.

Auto Start

This stirrer includes an optional Auto Start feature. When set, the stirrer will resume mixing after a power failure, when power has been removed by control equipment (i.e., PLC) or when the power has been cut (i.e., stirrer unplugged). When power is restored, the stirrer will automatically ramp up to the last set speed and resume mixing. The stirrer ships from the factory with Auto Start disabled.

To set the Auto Start feature first ensure the stirrer is powered off. Press and hold the Increase Button then press the Power Button. When active, Auto Start is indicated by:


BDC2002 - A solid LED dot in the bottom right corner of the display

BDC2010 - Flashing LED dots along the bottom of the display


To disable Auto Start first ensure stirrer is powered off. Press and hold the Decrease Button then press the Power Button.

Torque Capabilities


BDC2002 Torque Curve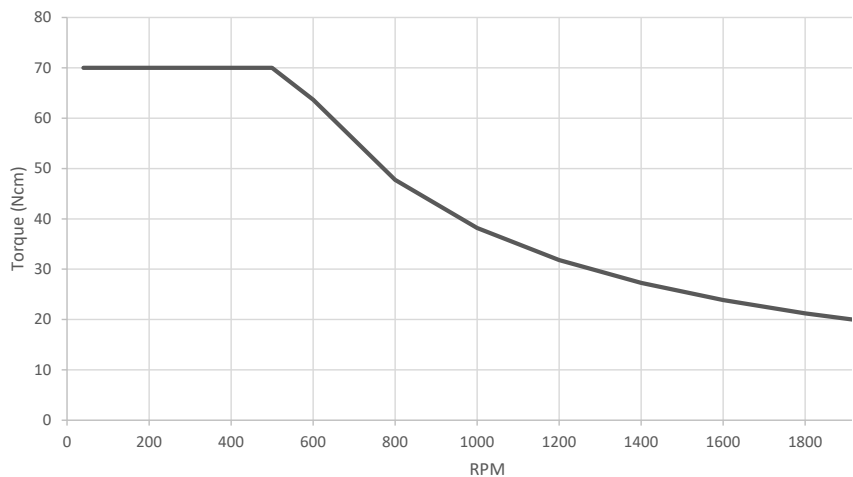

BDC2010 Torque Curve

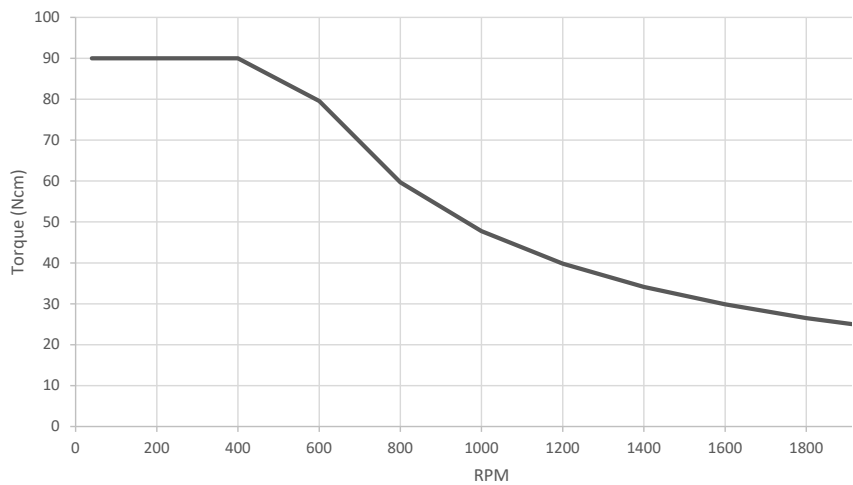
BDC2002 Dimensions

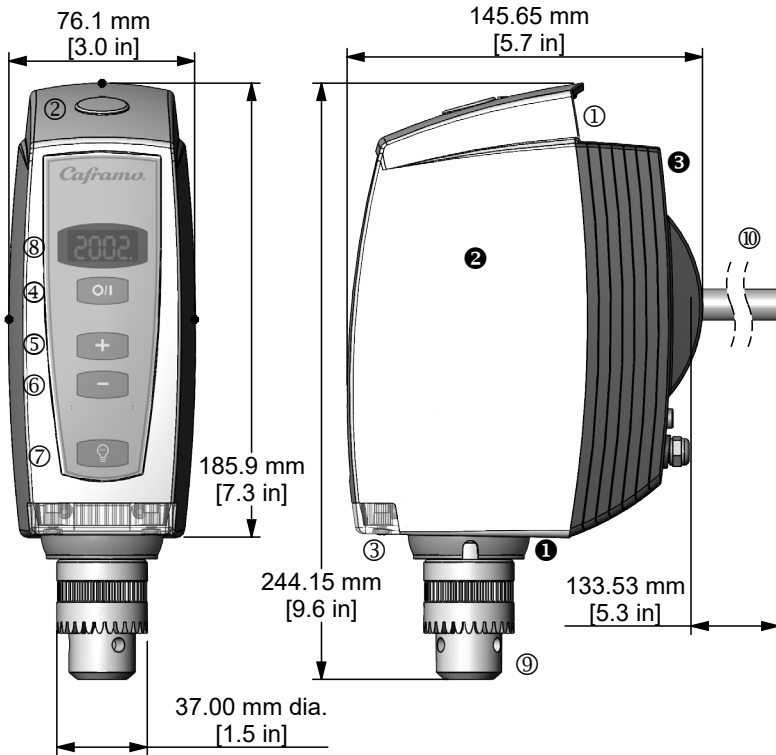- ① Ergonomic Carry Handle
- ② Rubber Gasket
- ③ Stirlight™
- ④ Button - Power on/off
- ⑤ Button - Increase rpm
- ⑥ Button - Decrease rpm
- ⑦ Button - Light on/off

- ⑧ Four Digit LED Display
- ⑨ Chuck
- ⑩ 1/2 in dia Support Rod
- ① Splash Guard
- ② Housing - Molded Front
- ③ Housing - Cast Back

BDC2010 Dimensions

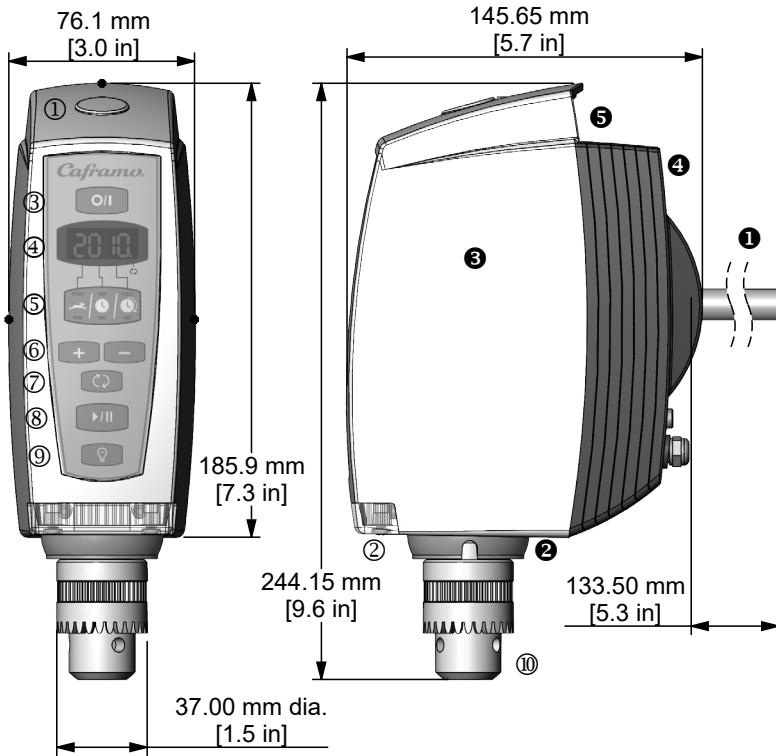- ① Rubber Gasket
- ② Stirlight™
- ③ Button - Power on/off
- ④ Four Digit LED Display
- ⑤ Multi-Function Button
- ⑥ Button - Increase/Decrease rpm
- ⑦ Button - Reverse direction
- ⑧ Button - Start/Pause
- ⑨ Button - Light on/off
- ⑩ Chuck
- ① 1/2 in dia Support Rod
- ② Splash Guard
- ③ Housing - Molded Front
- ④ Housing - Cast Back
- ⑤ Ergonomic Carry Handle

BDC2002 Technical Data

Model Number	BDC2002
Speed Range	40 - 2002 rpm
Speed Readout Accuracy	± 1 rpm or 3% of reading, 4 digit
Maximum Torque	70 Ncm, (100 in-oz)
Input To Power Supply	AC 100-240 volts, 50/60 Hz, 2.0 amps (Must use Caframo supplied power supply.)
Electrical Output from Power Supply to Stirrer	DC 24 volt, 3.1 amps
Mechanical Output Power	1/18 hp, 40 watts, (Brushless DC Motor)
Weight - Stirrer Only	3.0 kg (7 lbs)
Weight - Power Supply	0.5 kg (1 lbs)
Maximum Volume	25 litres (6.6 US gallons)
Maximum Viscosity	15,000 cps - e.g. sour cream
Molded Front Housing	Flame retardant, chemical resistant, high impact nylon. Recyclable (PA)..
Cast Back Housing	Epoxy coated zinc
Combined Housing	Splash Proof IP42
Keypad	Waterproof, chemical resistant polyester.
Stirlight™	5 Ultra Bright LEDs, 100,000 hour life.
Power Cord Length	183 cm (6 ft) from stirrer to power supply-DC. 183 cm (6 ft) from power supply to wall-AC.
Adjustable Chuck	Fits onto 13 mm (33/64 in) dia output shaft of stirrer. Adjusts up to 10 mm (3/8 in).

BDC2010 Technical Data

Model Number	BDC2010
Speed Range	40 - 2010 rpm
Functions	Countdown Timer Clockwise Counter Clockwise xRx Agitation
Speed Readout Accuracy	± 1 rpm or 3% of reading, 4 digit
Maximum Torque	90 Ncm, (127 in-oz)
Input To Power Supply	AC 100-240 volts, 50/60 Hz, 2.5 amps (Must use Caframo supplied power supply.)
Electrical Output from Power Supply to Stirrer	DC 24 volts, 3.75 amps
Mechanical Output Power	1/15 hp, 50 watts, (Brushless DC Motor)
Weight - Stirrer Only	3.0 kg (7 lbs)
Weight - Power Supply	0.6 kg (1.2 lbs)
Maximum Volume	25 litres (6.6 US gallons)
Maximum Viscosity	20,000 cps - e.g. chocolate syrup
Molded Front Housing	Flame retardant, chemical resistant, high impact nylon. Recyclable (PA).
Cast Back Housing	Epoxy coated zinc
Combined Housing	Splash Proof IP42
Keypad	Waterproof, chemical resistant polyester.
Stirlight™	5 Ultra Bright LEDs, 100,000 hour life.
Power Cord Length	183 cm (6 ft) from stirrer to power supply-DC. 183 cm (6 ft) from power supply to wall-AC.
Adjustable Chuck	Fits onto 13 mm (33/64 in) dia output shaft of stirrer. Adjusts up to 10 mm (3/8 in).

BDC2002 General Operating Instructions

Make sure the stirrer is securely mounted as specified in the installation instructions. Ensure the mixing paddle or impeller is securely attached at the desired position.

The stirrer is now ready for use to mix liquids and liquid/solid solutions. Appropriate impeller and container specifications must be determined by the operator, ensuring that all safety instructions are followed.

Keypad Functions

Power Button - Turns the stirrer motor on and off.

Increase Button - Increases the speed by one rpm. If pressed within five seconds of the stirrer being turned on it will stir at the last speed before powered off.

Decrease Button - Decreases the speed by one rpm. If pressed within five seconds of the stirrer being turned on it will stir at 40 rpm.

Light Power Button - Turns the light on and off. The stirrer must be plugged in for the light to function.

Safety is built into this stirrer. When it is turned on it pauses for 5 seconds before it starts to mix. The last speed used is displayed on the LED. The user has three options.

1. Do nothing. The stirrer will start to stir at 40 rpm after five seconds.
2. Press the increase button and the stirrer will start to stir at the last speed set before powered off. This speed is displayed.
3. Press the decrease button and the stirrer will start to stir at 40 rpm.

LED Display

Numeric value reflects speed (rpm). A solid LED dot indicates the Auto Start feature is enabled (see Auto Start). A “soft start” has been incorporated into the stirrer. When the increase or decrease buttons are pressed and held, the speed ramps smoothly so that getting to the speed you require is quick and easy.

Display Codes

OL Indicates an overload condition. The load is too great for the stirrer and must be reduced. To reset stirrer turn power off and back on.

OH Indicates an overheat condition. The operating conditions have caused the stirrer to overheat. Reduce stirrer loads and / or environmental temperature

to continue mixing. To reset stirrer turn power off, allow stirrer to cool and then turn back on.

BDC2010 General Operating Instructions

Make sure the stirrer is securely mounted as specified in the installation instructions. Ensure the mixing paddle or impeller is securely attached at the desired position.

The stirrer is now ready for use to mix liquids and liquid/solid solutions. Appropriate impeller and container specifications must be determined by the operator, ensuring that all safety instructions are followed.

Keypad Functions

Power Button - Turns the stirrer on and off.

Multi-Function Button - Sets three parameters - speed (rpm), countdown timer (minutes) and xRx agitation (seconds).

Increase Button - Increases the speed by one rpm or the countdown timer by one minute or the xRx agitation by one second.

Decrease Button - Decreases the speed by one rpm or countdown timer by one minute or the xRx agitation by one second.

Reverse - Reverses the direction that the impeller is turning. The impeller turns clockwise by default.

Start/Pause Button - Allows the stirrer to be paused and started again. Use this button to start the stirrer when ready.

Light Power Button - Turns the light on and off. The stirrer must be plugged in for the light to function.

LED Display - Numeric values reflect speed (rpm), countdown timer (minutes) or xRx agitation (seconds). Solid LED dots indicate which function is selected. Flashing LED dots indicate the Auto Start feature is active (see Auto Start). The LED dots do not act as numeric decimal points.

① LED dot (solid) when speed is displayed.

② LED dot (solid) when time is displayed.

③ LED dot (solid) when cycle time is displayed.

↻ LED dot (solid) when stirrer is turning counter clockwise.

Clockwise Mixing

Turn the stirrer on by pressing the power button. The display will read 40 rpm (or last set speed when Auto Start is enabled) and the speed LED dot will be illuminated. Increase to the desired speed and press the start/pause button. The stirrer will start turning in a clockwise direction. The speed LED dot will be illuminated (solid) and the speed value in rpm will be displayed.

Counter Clockwise Mixing

Set by following the clockwise stirring instructions above. Press the reverse button. The reverse button can be pressed while the stirrer is running to mix in a counter clockwise direction. The LED dot above the reversing icon will be illuminated (solid) when mixing in a counter clockwise direction.

Countdown Timer

Turn the stirrer on by pressing the power button. Set the desired speed by pressing the “+” button. Press the multifunction button until the LED dot for time is illuminated. Press the “+” to set the countdown timer in minutes. Press the start/pause button to start mixing. Press the multifunction button to toggle the display between speed, time and cycle. Stirring stops when time is completed. Three beeps will sound. Note: Countdown Timer value cannot be changed during operation.

xRx Agitation Cycle

Turn the stirrer on by pressing the power button. Set speed as above and countdown timer if desired. Press the multifunction button once more until the indicator light for cycle is illuminated. Press the “+” to set the number of seconds the impeller will rotate in one direction before changing directions. Press the start/pause button. Note: xRx Agitation Cycle value cannot be changed during operation.

For example: If set for 5 seconds the impeller will turn clockwise for 5 seconds at the set speed then stop and change to turn counter clockwise for 5 seconds. This pattern will continue until the pause button is pressed or until the time is completed (if set).

The display will show the cycle seconds counting down during each direction unless speed or the countdown timer is selected.

When the stirrer is powered on, the default speed is 40 rpm, countdown timer is 0 minutes, and the xRx agitation cycle is 0 seconds. Previously set values will be displayed if Auto Start is enabled.

The stirrer will remember the original values set for speed, the countdown timer and the xRx agitation cycle so that when the start/pause button is pressed again the operations will repeat.

Speed can be changed (increased or decreased) while the stirrer is running. Press

the multifunction button so the speed is displayed. Increase or decrease speed by pressing the “+” or “-” buttons.

The countdown timer and the xRx agitation cycle can be changed only when the stirrer is paused and the indicator lights are solid.

The multifunction button can be pressed when the stirrer is running to view the status of each function.

For safety, a “soft start” has been incorporated into the stirrer. The stirrer will ramp up to reach the set speed to minimize splashing.

Display Codes

OL Indicates an overload condition. The load is too great for the stirrer and must be reduced. Press the power button, enter in new parameters and start.

OH Indicates an overheat condition. The operating conditions have caused the stirrer to overheat. Reduce stirrer loads and/or environmental temperature to continue mixing. To reset stirrer turn power off, allow stirrer to cool and then turn back on.

Cleaning and Service

Cleaning

Unplug stirrer prior to cleaning. Using a mild soap and water solution, wipe the housing and control panel of the stirrer with a damp cloth. Do not use chlorine bleach, chlorine-based cleanser, abrasives, ammonia, steel wool, or scouring pads with metal content or similar harsh solvents or abrasives. These may damage the surface of the stirrer.

Service

Do not attempt to service or repair stirrer. If the stirrer housing is opened, the warranty becomes void. Contact Supplier for return authorization and return instructions. Clean the stirrer so that it is free from any materials which may be harmful. Provide an SDS or other safety data sheets for all chemicals used in conjunction with this stirrer where appropriate. Place the stirrer and its parts into the original packing to protect and prevent damage during transport. Follow shipping directions provided by Caframo when return is authorized and RMA issued.

CAUTION:

 Failure to clean, service, and handle the stirrer as outlined above can cause it to fail which could impair the safety of the user.

Warranty and Product Liability

Warranty

Caframo warrants the herein described product of its manufacture for three years from receipt of the equipment by the purchaser, against defects in material and workmanship. Warranty registration is available on the Service and Support page of our website (CaframoLabSolutions.com).

This limited warranty covers parts and labour. If the stirrer housing is opened, the warranty becomes void.

This warranty may only be altered by a specifically published amendment. No individual has authorization to alter the provisions of this warranty policy or its amendments. Expendable items are not covered by this warranty. Damages due to corrosion or accidental breakage are not covered.

Product Liability

Under no circumstances shall Caframo be liable for indirect, consequential, or special damages of any kind.

Caframo reserves the right to make technical changes without prior notice.

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