

PolySep™

Polymer-Based Aqueous GFC Columns for the Separation of Polymers, Oligomers, and Oligosaccharides

Extremely cost-effective

High efficiencies

Good mechanical strength

 **phenomenex**[®]
...breaking with traditionSM




Dextran

Column: PolySep-GFC-P4000
Dimensions: 300 x 7.8 mm
Part No.: 00H-3144-K0
Mobile Phase: Water
Flow Rate: 0.8 mL/min
Detection: RI

Sample: 20 µL injected

1. Dextran	500	kDa
2. Dextran	110	kDa
3. Dextran	70	kDa
4. Dextran	40	kDa
5. Dextran	15	kDa
6. Dextran	6	kDa
7. Ethylene Glycol	62.07	Da

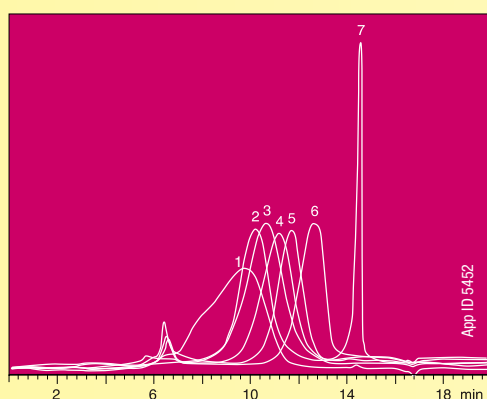


Polyethylene Standards

Column: PolySep-GFC-P4000
Dimensions: 300 x 7.8 mm
Part No.: 00H-3144-K0
Mobile Phase: Water
Flow Rate: 0.8 mL/min
Detection: RI

Sample: 20 µL injected

1. Polyethylene Oxide	305.4	kDa
2. Polyethylene Oxide	95.7	kDa
3. Polyethylene Oxide	31.8	kDa
4. Polyethylene Glycol	19.7	kDa
5. Polyethylene Glycol	3.4	kDa
6. Ethylene Glycol	62.07	Da

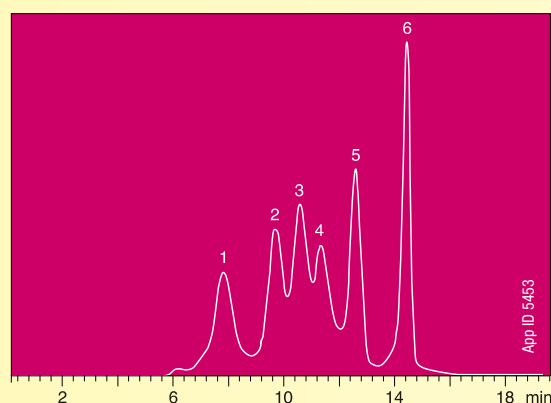


Polyethylene Oxide/Polyethylene Glycol

Column: PolySep-GFC-P3000
Dimensions: 300 x 7.8 mm
Part No.: 00H-3143-K0
Mobile Phase: Water
Flow Rate: 0.8 mL/min
Detection: RI

Sample: 10 µL injected

1. Polyethylene Oxide	62.6	kDa
2. Polyethylene Oxide	31.1	kDa
3. Polyethylene Glycol	19.7	kDa
4. Polyethylene Glycol	6.8	kDa
5. Polyethylene Glycol	3.4	kDa
6. Polyethylene Glycol	1.5	kDa
7. Ethylene Glycol	62.07	Da

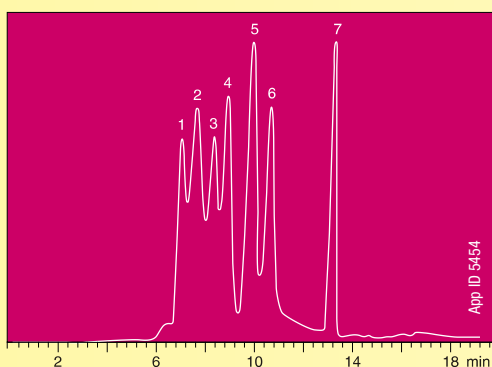

Aqueous SEC 2 Standard (Part No. AL0-3043)


Column: PolySep-GFC-P1000
Dimensions: 300 x 7.8 mm
Part No.: 00H-3141-K0
Mobile Phase: Water
Flow Rate: 0.8 mL/min
Detection: RI

Sample: 3 µL injected

1. Ethylene Glycol

Analyst Note: This standard is used to test the following PolySep columns: GFC-P 1000, 2000, 3000, 4000, 5000, 6000, and Linear, as well as other aqueous-soluble GFC columns.

NOTE: Separation conditions and results may vary for your column. Contact your Technical Representative.

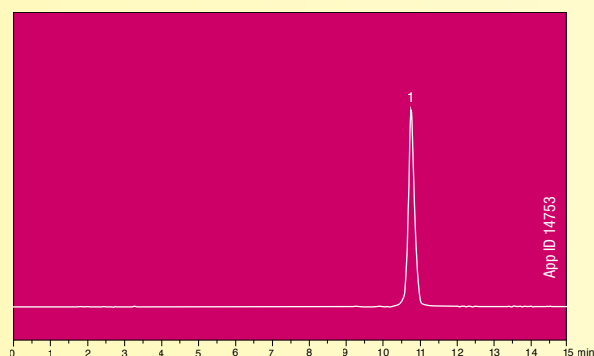

Polymer-Based Aqueous GFC Columns for the Separation of Polymers, Oligomers, and Oligosaccharides

The PolySep material undergoes rigorous quality control tests, from the initial stages of testing of the starting monomers to the final product. There are at least 25 steps of quality assurance during the entire procedure. The packed column then undergoes at least five additional tests, including a batch test for the manufactured materials. Each column is then tested for column efficiency and peak symmetry and ships with a QC chromatogram. This ensures long-lasting columns with very high efficiencies.

PolySep-GFC-P Technical Data and Specifications

Phase:	1000	2000	3000	4000	5000	6000	Linear
Exclusion Limits in Daltons:							
PEG	2 x 10 ³	9 x 10 ³	5 x 10 ⁴	2 x 10 ⁵	2 x 10 ⁶	1 x 10 ⁷	1 x 10 ⁷
Pullulans	3.5 x 10 ³	1 x 10 ⁴	1 x 10 ⁵	3.5 x 10 ⁵	4 x 10 ⁶	2 x 10 ⁷	2 x 10 ⁷
Separation Range (Da)	20 - 3K	100 - 10K	250 - 75K	3K - 400K	50K - 2M	100K - 15M	1K - 10M
Typical Efficiency Plates/meter	22,000	50,000	32,000	32,000	32,000	32,000	32,000
Maximum Organic Modifier:							
Methanol	20 %	95 %	70 %	70 %	70 %	70 %	70 %
Acetonitrile	20 %	70 %	70 %	70 %	70 %	70 %	70 %
pH Range	3.0 to 12.0						
Maximum Flow Rate	Depends on backpressure, do not exceed 1000 psi						
Column Hardware	Stainless steel or PEEK (Biocompatible hardware available upon request)						
Temperature	4 to 60 °C						
Maximum Salt	Maximum allowed 0.5 M with a flow rate not to exceed 0.5 mL/min						
Storage	For overnight, pump water at 0.2 mL/min, for longer storage use 0.05 % NaN ₃ in water or 10 % methanol in water						
General	A guard column is recommended to improve column life						

Ordering Information

Stainless Steel Columns (mm)

	Analytical	Guards
Phases	300 x 7.8	35 x 7.8
1000	CH0-9226	CH0-9225
2000	CH0-9227	CH0-9225
3000	CH0-9228	CH0-9225
4000	CH0-9229	CH0-9225
5000	CH0-9230	CH0-9225
6000	CH0-9231	CH0-9225
Linear	CH0-9232	CH0-9225

App ID 14753

Aqueous SEC 2

(For PolySep GFC-P and other aqueous-soluble analysis columns)

Part No.: AL0-3043

Price:

Unit quantity: 2 mL

Contains: Ethylene Glycol

Diluent: Water

Test Conditions

Mobile Phase: Water

Flow Rate: 0.8 mL/min

Injection Volume: 15 µL

Detection: RI

Economical alternative to

- Tosoh Bioscience® TSKgel® PW & PWxl
- Agilent® PL aquagel-OH™
- Waters® Ultrahydrogel™
- Shodex® OHpak SB-800 HQ series
- PSS SUPREMA®



Polymer-Based Aqueous GFC Columns for the Separation of Polymers, Oligomers, and Oligosaccharides

Australia

t: +61 (0)2-9428-6444
f: +61 (0)2-9428-6445
auinfo@phenomenex.com

Austria

t: +43 (0)1-319-1301
f: +43 (0)1-319-1300
anfrage@phenomenex.com

Belgium

t: +32 (0)2 503 4015 (French)
t: +32 (0)2 511 8666 (Dutch)
f: +31 (0)30-2383749
beinfo@phenomenex.com

Canada

t: +1 (800) 543-3681
f: +1 (310) 328-7768
info@phenomenex.com

China

t: +86 (0)20 2282-6668
f: +86 (0)20 2809-8130
chinainfo@phenomenex.com

Denmark

t: +45 4824 8048
f: +45 4810 6265
nordicinfo@phenomenex.com

Finland

t: +358 (0)9 4789 0063
f: +45 4810 6265
nordicinfo@phenomenex.com

France

t: +33 (0)1 30 09 21 10
f: +33 (0)1 30 09 21 11
franceinfo@phenomenex.com

Germany

t: +49 (0)6021-58830-0
f: +49 (0)6021-58830-11
anfrage@phenomenex.com

India

t: +91 (0)40-3012 2400
f: +91 (0)40-3012 2411
indiainfo@phenomenex.com

Ireland

t: +353 (0)1 247 5405
f: +44 1625-501796
eireinfo@phenomenex.com

Italy

t: +39 051 6327511
f: +39 051 6327555
italiainfo@phenomenex.com

Luxembourg

t: +31 (0)30-2418700
f: +31 (0)30-2383749
nlinfo@phenomenex.com

Mexico

t: 001-800-844-5226
f: 001-310-328-7768
tecnicomx@phenomenex.com

The Netherlands

t: +31 (0)30-2418700
f: +31 (0)30-2383749
nlinfo@phenomenex.com

New Zealand

t: +64 (0)9-4780951
f: +64 (0)9-4780952
nzinfo@phenomenex.com

Norway

t: +47 810 02 005
f: +45 4810 6265
nordicinfo@phenomenex.com

Puerto Rico

t: +1 (800) 541-HPLC
f: +1 (310) 328-7768
info@phenomenex.com

Spain

t: +34 91-413-8613
f: +34 91-413-2290
espinfo@phenomenex.com

Sweden

t: +46 (0)8 611 6950
f: +45 4810 6265
nordicinfo@phenomenex.com

United Kingdom

t: +44 (0)1625-501367
f: +44 (0)1625-501796
ukinfo@phenomenex.com

USA

t: +1 (310) 212-0555
f: +1 (310) 328-7768
info@phenomenex.com

All other countries Corporate Office USA

t: +1 (310) 212-0555
f: +1 (310) 328-7768
info@phenomenex.com



www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department at international@phenomenex.com

Terms and Conditions

Subject to Phenomenex Standard Terms & Conditions, which may be viewed at www.phenomenex.com/TermsAndConditions.

Trademarks

Phenomenex is a registered trademark, PolySep is a trademark of Phenomenex. Waters is a registered trademark, Ultrahydrogel is a trademark of Waters Corporation. Tosoh Bioscience and TSKgel are registered trademarks of Tosoh Corporation. Agilent is a registered trademark, PL aquagel-OH is a trademark of Agilent Technologies, Inc. Shodex is a registered trademark of Showa Denko K.K. Suprema is a registered trademark of PSS.

Disclaimer

Phenomenex is not affiliated with Waters Corporation, Tosoh Corporation, Agilent Technologies, Inc, Showa Denko K.K or PSS.